

RECOMMENDED DILUTIONS

Kit name	Kit code	Juice	Finished wine
Acetic acid	4A100	1 in 5	1 in 5
	4A105		
PAAN	4A110	1 in 2	1 in 2
Ammonia	4A120	1 in 10	1 in 5
Citric acid	4A126	1 in 2	
Gluconic acid	4A130	1 in 10	
Glucose and fructose	4A140	1 in 50	1 in 2
	4A145		
Malic acid	4A160	1 in 5	1 in 2
	4A165		
Sucrose	4A180		
free SO ₂	4A199		
Total SO ₂	4A200		

HOW TO MAKE 1 mL DILUTED SAMPLE

Dilution needed	Volume of sample μ L	Volume of water (μ L)
1 in 2	500	500
1 in 5	200	800
1 in 10	100	900
1 in 50	20	980

Disclaimer – These values are indicative values based on results obtained from a NATA accredited laboratory with calibrated equipment.